Bysameyee USB Digital Microscope User Manual

Email Support: bysameyee@outlook.com

Web: <https://www.bysameyee.com/microscope>

- Thank you for purchasing our USB Digital Microscope and taking the time to read this guideline. We recommend you reading this manual **FIRST** to get the best out of this unit.
- Operation system: Windows 7 8 10/Vista/XP/2000, Mac OS X 10.5 or above, Android mobiles (with OTG&UVC function), Linux, Chrome OS.
- **Note:** The Zoom button only works on the Windows system with Amcap software.

Important!

Remove the lens cover before using the microscope.

1. Windows

The microscope is plug and play for Windows10/11 users, you can just plug in the microscope and run the system built-in webcam to use it. (The zoom and snap button won't work with win10/11 webcam)

All Windows users(win7/8/10/11) can download the Amcap software online: www.bysameyee.com/microscope


 [Amcap.zip\(3.34MB\) | Windows XP, Vista, 7, 8, 10 | 32/64 bit](#)

After downloaded, right-click the amcap.zip to extract it to your PC.


Then open the Amcap folder and double-click the setup.exe to install.

Name	Date modified	Type	Size
 data2.cab	3/18/2017 7:33 PM	Cabinet File	1 KB
 engine32.cab	3/18/2017 7:33 PM	Cabinet File	450 KB
 glavcam.cat	3/20/2017 4:26 PM	Security Catalog	12 KB
 glavcam.cer	3/18/2017 7:33 PM	Security Certificate	1 KB
 glavcam.inf	3/20/2017 4:24 PM	Setup Information	16 KB
 glavcam.sys	3/18/2017 7:33 PM	System file	2,713 KB
 glprop.ax	3/18/2017 7:33 PM	AX File	96 KB
 layout.bin	3/18/2017 7:33 PM	BIN File	1 KB
 monitorpad.exe	3/18/2017 7:33 PM	Application	68 KB
 monitorpad.ini	3/18/2017 7:33 PM	Configuration sett...	1 KB
 ProductName.ini	3/19/2017 8:52 AM	Configuration sett...	1 KB
 readme.txt	3/19/2017 8:54 AM	Text Document	3 KB
 setup.exe	3/18/2017 7:33 PM	Application	114 KB
 setup.ibt	3/18/2017 7:33 PM	IBT File	355 KB
 setup.ini	3/18/2017 7:33 PM	Configuration sett...	1 KB
 setup.inx	3/18/2017 7:33 PM	INX File	204 KB

Click “Next –Next -Install –Finish” to install the driver. And then select “No, I will restart my computer later”. Plug the microscope into the computer and open the Amcap driver. The default camera may not be the USB microscope, then please select GL USD2.0 UVC Camera Device from the “Devices” drop down menu. If you can't find this device, contact us and we will assist you to update it.

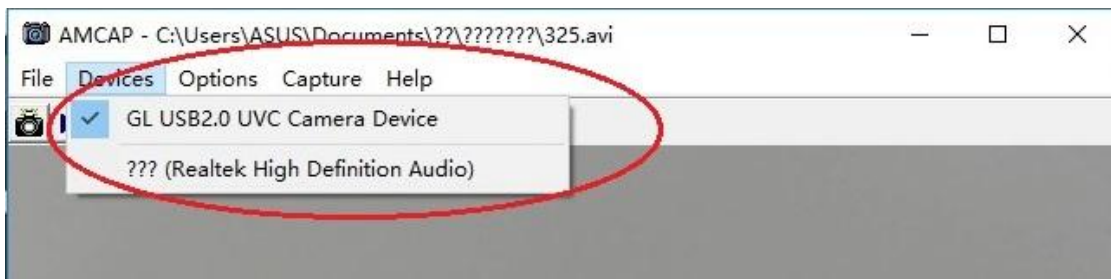
Note: If the Amcap software occurs error “Sorry, you have no video capture hardware”, please close the Amcap and right-click the Amcap.exe and select Properties – Compatibility and select “Run this program as an administrator”.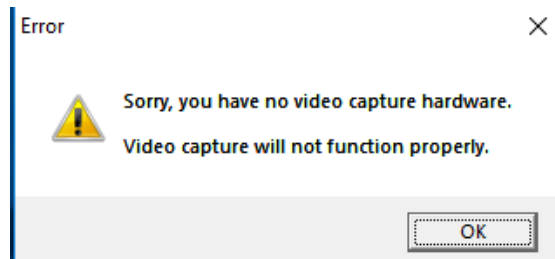

You need to set up the snapshot location for pictures and videos on Amcap. Pictures: Select “Capture” – Snapshot Location – Find or new a folder and click “Open”.

Videos: Click “File” – “Set Capture File”. Find and select an Avi format video file. Click “Open”.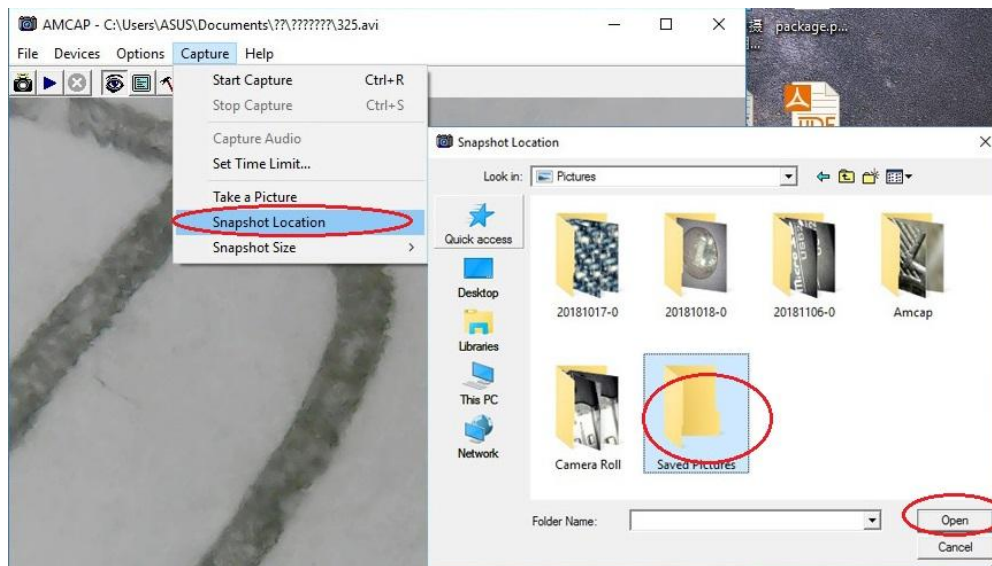

Measurement Tool Download (Optional):


Visit www.bysameyee.com/microscope and download:

When it is completed, right-click the Coolingtech.zip file and select Extract All. Open the CoolingTech folder and double click the CoolingTech.exe to use the measurement software. Click Run if it occurs security warning.

Note: The SNAP and Zoom button can't be used in the measurement software.


Note: Win10 users please set up the privacy setting on computer to let apps access the camera.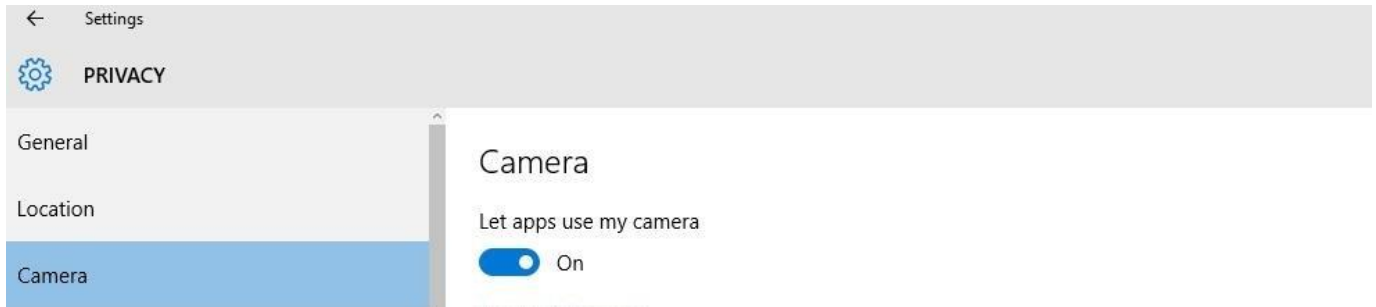

Get the newest update and alternatively applications on our website www.bysameyee.com/microscope

2. Mac

Mac users can use the built-in camera app “Photo Booth” to observe and capture pictures or videos. (Images are mirrored left-to-right)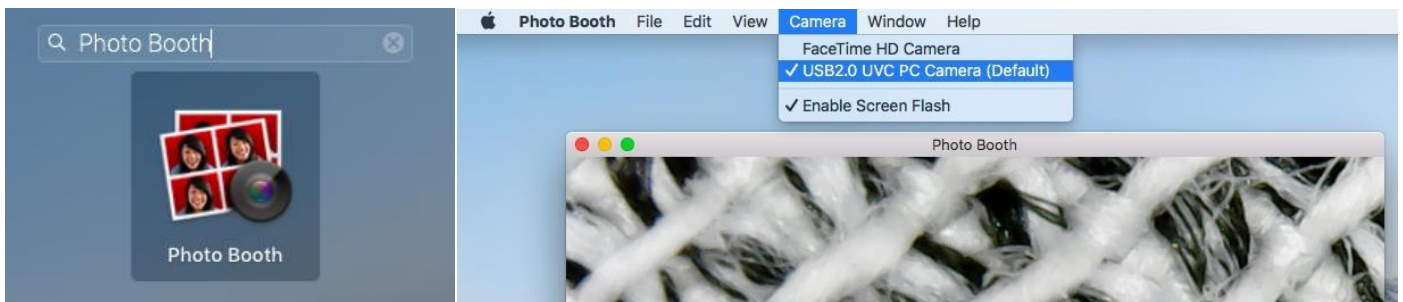


You can also install the backup camera app on www.bysameyee.com/microscope

Select one link to download the driver according to your Mac version:

↓ [Digital_Viewer_3.3.30.pkg\(8.66MB\) | macOS \(10.13.4 and later\) | 64 bit](#)

↓ [Digital_Viewer_3.1.06.dmg\(2.67MB\) | Older macOS / OS X versions \(10.5 and later\) | 32 bit](#)

Plug the microscope into the Mac computer and run the Digital Viewer app, click “Settings” – “Device” to select “USB2.0 UVC PC Camera” on the dropdown list, then click "Apply" to make the microscope work.

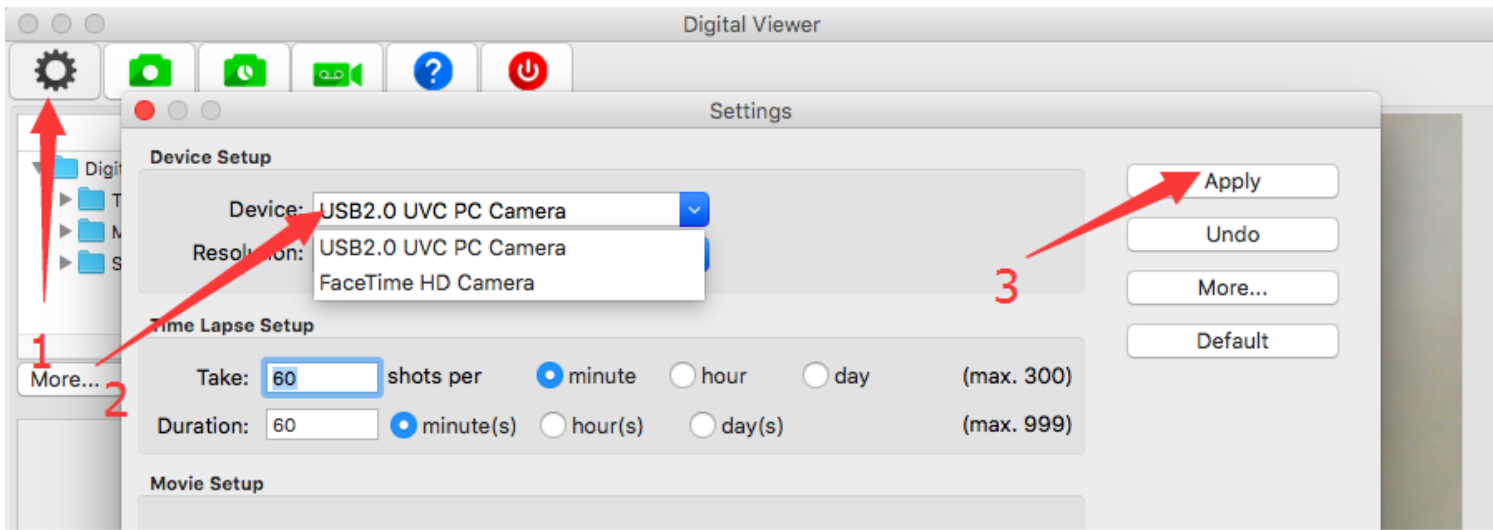3. Android

Before using, please make sure your Android phone supports **OTG and UVC function** and activated your OTG function (If optional).

Micro-USB phone users please flip a piece from the USB plug.

Type-C phone users please connect the Type-C adapter to the USB plug.

Apps Download:

Please download one of these apps on Google Play Store.

USB Camera, Otg View, AN98, xploview

Please note that all the apps we offer are totally **FREE**. Please contact us bysameyee@outlook.com when it requires paying for it.

Errors:

1) The microscope light doesn't light up when it's plugged in.

2) The light is on, but it shows a black screen, no picture.

Solution: Change USB adapter and plug-in again; Reinstall App or try to install another App.

Note: The Zoom button doesn't work on the phone app.

4. Chrome OS

This microscope can be used with the default Camera app that comes with Chrome. Some Chrome machines with built-in cameras may need to have the device pointer changed. This can be done by going to **Settings > Advanced Settings > Privacy > Site Settings > Camera**, and change the device in the dropdown to "USB Microscope". Open the Camera app on your Chromebook. You'll find it under the launcher menu. Tap the "Search" button on the keyboard and search for "Camera". Alternatively, click the "All Apps" button and look for the camera icon.

Get the newest update on our website www.bysameyee.com/microscope

5. Linux

Raspberry Pi: Working with the current version of Raspbian Jessie.

We have tested the microscope with Linux utilizing GTK+ UVC Viewer by using the following terminal commands:

“sudo apt-get install guvcview” “guvcview”

6. iPhone/iPad(Not support)

Since Apple Company blocks the data transfer function, this microscope cannot work with iPhone/iPad. If you ordered it by mistake, we can offer a special discount for you to order a wireless iPhone microscope. Please contact bysameyee@outlook.com

Q&A


Q: When running the Amcap driver, it occurs “Error: Cannot create video capture filter”.

A: Right-click the Amcap.exe and select **Properties – Compatibility** and select **run this program as an administrator**. Click **“OK”**.

Q: The software requires me to pay for it?

A: Don’t install it. Please contact us for a free version.

Q: Occur error as below show, “Access to the path ‘C:\Program Files (x86)\CoolingTech\Capture’ is denied when running the Coolingtech application.

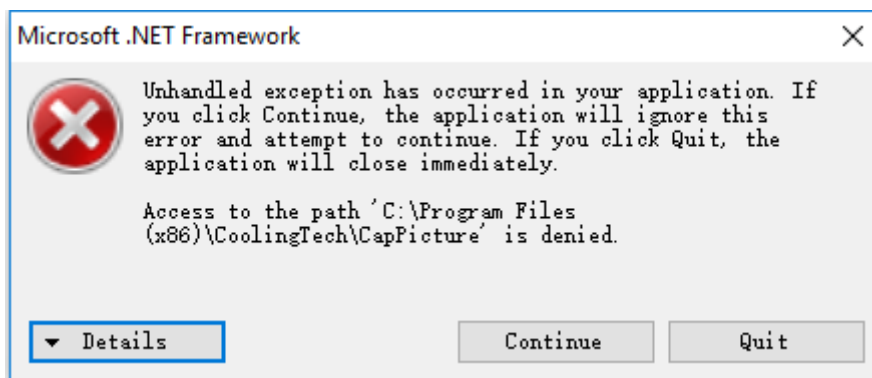A: Create 2 new folders and rename them as **“CapPicture”** and **“CapVideo”**.

Copy these 2 folders and paste them to **C:\Program Files (x86)\CoolingTech**.

If you install the Coolingtech in another path, just follow the reminder message to find the path.

Name	Date modified	Type	Size
CapPicture	9/14/2018 3:43 PM	File folder	
CapVideo	9/14/2018 3:43 PM	File folder	
CoolingTech	3/16/2017 12:25 AM	Application	2,509 KB
CoolingTech	3/16/2017 12:25 AM	ICO File	14 KB
DirectShowLib-2005.DLL	3/16/2017 12:25 AM	Application extens...	296 KB
Interop.Outlook.DLL	2/24/2017 8:04 PM	Application extens...	392 KB

Q: Signature issue.(Code 52)

A: Please plug the microscope into the computer and right-click “[This PC](#)” – “[Properties](#)” and select “[Device Manager](#)” to open the device manager. Find the “[GL USB2.0 UVC Camera Device](#)” on the Imaging device or Cameras menu. Right-click and select “[Uninstall device](#)” and select “[Delete the driver software for this device](#)”. Click “[Uninstall](#)” “[Yes](#)” to restart the computer. Open the Amcap to use the microscope once the computer is restarted.

Please visit <https://www.bysameyee.com/microscope> for newest Q&A for this microscope.

To save your time, don't hesitate to email us bysameyee@outlook.com if there's any issue with the order or even you require a **discount** from us.

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Name	Date modified	Type	Size
data2.cab	3/18/2017 7:33 PM	Cabinet File	1 KB
engine32.cab	3/18/2017 7:33 PM	Cabinet File	450 KB
glavcam.cat	3/20/2017 4:26 PM	Security Catalog	12 KB
glavcam.cer	3/18/2017 7:33 PM	Security Certificate	1 KB
glavcam.inf	3/20/2017 4:24 PM	Setup Information	16 KB
glavcam.sys	3/18/2017 7:33 PM	System file	2,713 KB
glprop.ax	3/18/2017 7:33 PM	AX File	96 KB
layout.bin	3/18/2017 7:33 PM	BIN File	1 KB
monitorpad.exe	3/18/2017 7:33 PM	Application	68 KB
monitorpad.ini	3/18/2017 7:33 PM	Configuration sett...	1 KB
ProductName.ini	3/19/2017 8:52 AM	Configuration sett...	1 KB
readme.txt	3/19/2017 8:54 AM	Text Document	3 KB
setup.exe	3/18/2017 7:33 PM	Application	114 KB
setup.ibt	3/18/2017 7:33 PM	IBT File	355 KB
setup.ini	3/18/2017 7:33 PM	Configuration sett...	1 KB
setup.inx	3/18/2017 7:33 PM	INX File	204 KB

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